Agilent SPS 4 and 6 Packing/Unpacking Guide

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Unpacking Instructions/ 开箱说明 / 開梱手順

Step	Action
------	--------

WARNING

Heavy

Danger to hands and feet. When unpacked the autosampler is heavy. To avoid injury to personnel or damage to equipment, always use at least two people to lift the autosampler into position.

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-
- | | |
|---|---|
| 1 | Before accepting delivery, check the shipping box for any external signs of damage. If the box shows any sign of damage, contact the carrier immediately. Any damage should be noted on the carrier's way-bill. |
|---|---|
-

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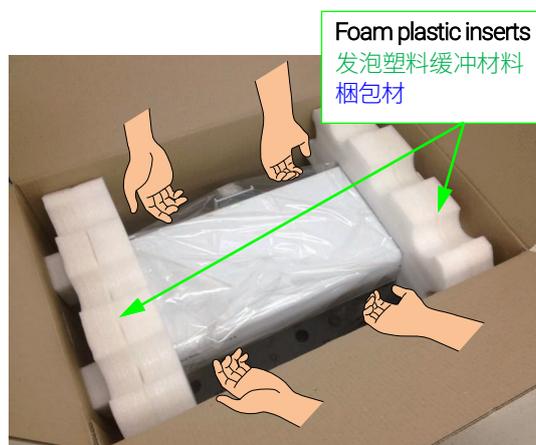
Step Action

- 2 Unpack the shipping box and lift the unit and all accessories from the box.
One person lifts from the front and one from the rear of the unit. Bend at the knees if lifting from a low position. Remove these foam plastic inserts before lifting unit.

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- 3 Place the unit on a suitable bench and check all of the items you have received.

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If you require detailed instructions on unpacking, refer to "Packing Instructions" on Page 3 and follow the steps in reverse order.

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NOTE

Keep the shipping box and packaging. If you need to return the autosampler during the warranty period, it is a requirement of the warranty that the equipment is repacked in the original shipping box.

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CAUTION

Inspect all items for damage. Any damage or missing items should be reported to Agilent immediately. Cite the part number and serial number of your autosampler.

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Packing Instructions/ 包装说明 / 梱包手順

If you need to return the autosampler to Agilent Technologies for repair, follow the steps below to repack the equipment.

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- | | |
|---|--|
| 1 | <ul style="list-style-type: none">• Move the sample probe mount all the way down.• Remove all accessories from the autosampler except for the spill tray. |
|---|--|

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- | | |
|---|---|
| 2 | <ul style="list-style-type: none">• Cover the probe arm with a suitable plastic bag.• Place the probe arm and plastic insert against the left end frame and wrap stretch wrap around the top half to secure it in place. Wrap the plastic insert to prevent it from moving vertically. |
|---|---|

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Step Action

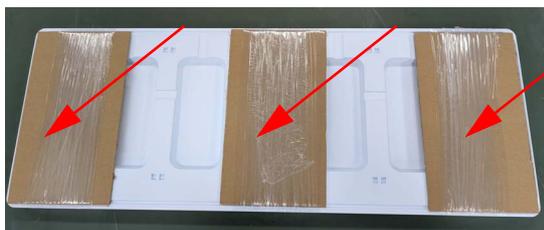
- 3
- For the SPS 4, to reinforce the spill tray, place cardboard in the center of the tray and secure it with stretch wrap.
 - For the SPS 6, to reinforce the spill tray, place cardboard in the center and sides of the tray and secure it with stretch wrap.
 - Turn over the spill tray.
 - Place any accessories on top of the spill tray and secure them in place by wrapping them with stretch wrap.

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
- 4
- Place the spill tray upside down on the top of the base bars.
 - Use plastic inserts and stretch wrap to secure the spill tray through the gaps in the end frames.
 - For the SPS 6, apply additional stretch wrap from the front to the rear of the spill tray to secure both the left and right sides of the accessories.

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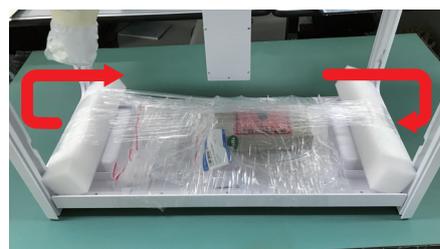Step Action

- 5 • Use stretch wrap to secure the Pressure Plates of the Peristaltic Pump so that they are not dislodged during shipping.

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- 6 • Place the unit inside the plastic bag and seal it with packing tape.

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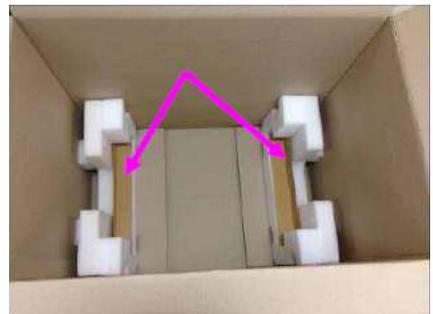
- 7 • Place the bottom plastic inserts at the bottom of the box along with a 400 mm piece of angled cardboard bracing.

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- 8 • Place the unit inside the box making sure the angled cardboard braces are still in place, by at least two persons.
• Place cardboard braces on top of the unit at both ends. Ensure there is no tape under the cardboard and that any markings on the cardboard are not facing the top of the unit.

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Step Action

- 9
- Place plastic inserts on top of the unit at both ends making sure the angled cardboard braces are still in place.
 - Add any relevant documentation in a sealable plastic bag taped to the top of the unit. Make sure the Unpacking Instructions are clearly visible when the box is opened.

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- 10
- For the SPS 4, close and seal the shipping box using packing tape and reinforced tape in an 'H' pattern.
 - For the SPS 6, close and seal the shipping box using packing tape and reinforced tape in a 'U' pattern.
 - Place labels as shown and make sure the orientation of the "This Way Up" labels are correct. Labels must show 'THIS WAY UP', 'FRAGILE', 'KEEP OUT OF RAIN', and 'HEAVY'.
 - Cut the tape over the handles and tuck back up inside the box.

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