

Agilent SPS 4 Autosampler

Cart Assembly Instructions

This document provides instructions on how to assemble the Agilent Autosampler cart.





Specifications

- Maximum of 120 kg (265 lbs) per shelf
- Maximum total loading capacity 250 kg (551 lbs)
- 100 mm diameter casters with two brakes and two swivels

CAUTION Only use the supplied rubber mallet to assemble the cart.


Do not apply excessive force when hammering the shelves onto the posts.

Parts List

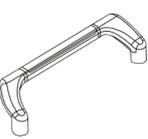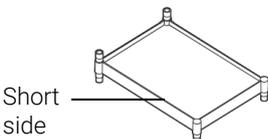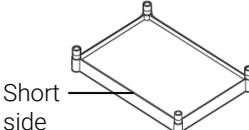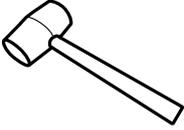- | | | | |
|---|---|--|---|
|  |  | 
Short side |  |
| 1 Handle – Qty 1 | 2 Separator – Qty 2 | 3 Top shelf – Qty 1 | 4 Washer – Qty 2 |
|  |  | 
Short side |  |
| 5 Screws – Qty 2 | 6 Posts – Qty 4 | 7 Bottom shelf – Qty 1 | 8 Split washer – Qty 4 |
|  |  |  |  |
| 9 Casters with lock – Qty 2 | 10 Swivel casters – Qty 2 | Wrench | #2 Phillips screwdriver |
| | | |  |
| | | | Rubber mallet |


Agilent SPS 4 Autosampler Cart Assembly Instructions

Step 1

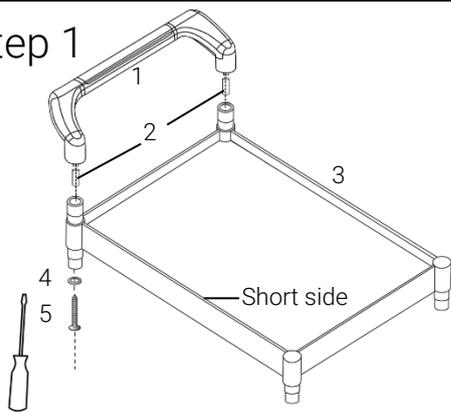


- a Insert separators (2) onto both handle (1) stand-offs.
- b Slide the handle into the top shelf (3).
- c Firmly secure both sides with a washer (4) and screw (5).

Step 2

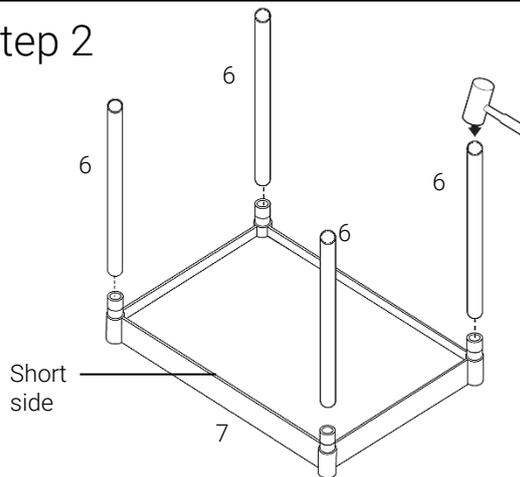


- a Push, and then hammer, the posts (6) onto the base (7).

Step 3

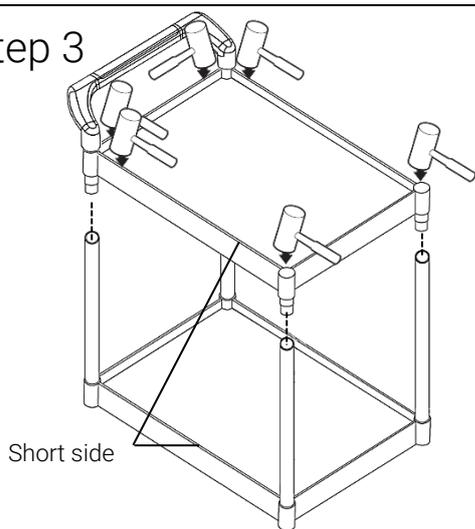

- a Push the top shelf into the posts and then evenly hammer each side until fully seated.


Step 4

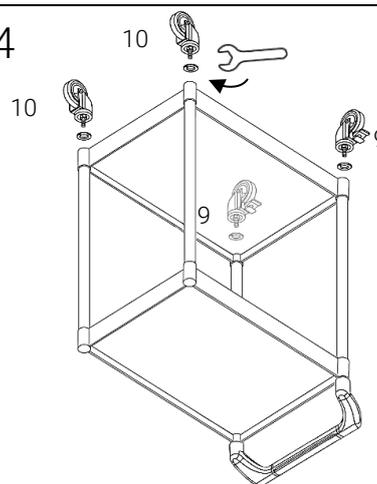

- a Flip the trolley upside down.
- b Screw in the washers (8) and casters (9 and 10). Use the wrench to tighten.

Step 5

- a Lock the casters and then carefully flip the cart right side up.
- b Unlock the casters to move the cart.

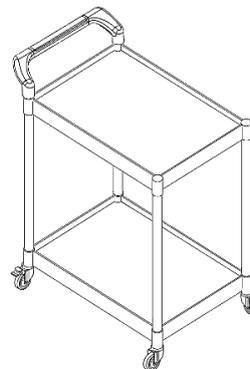

This information is subject to change without notice.

Part Number: G8410-90060

Edition 8/25
Issue 2
Printed in USA

© Agilent Technologies, Inc. 2025

Agilent Technologies Australia (M) Pty Ltd
679 Springvale Road
Mulgrave, VIC 3170, Australia



G8410-90060

